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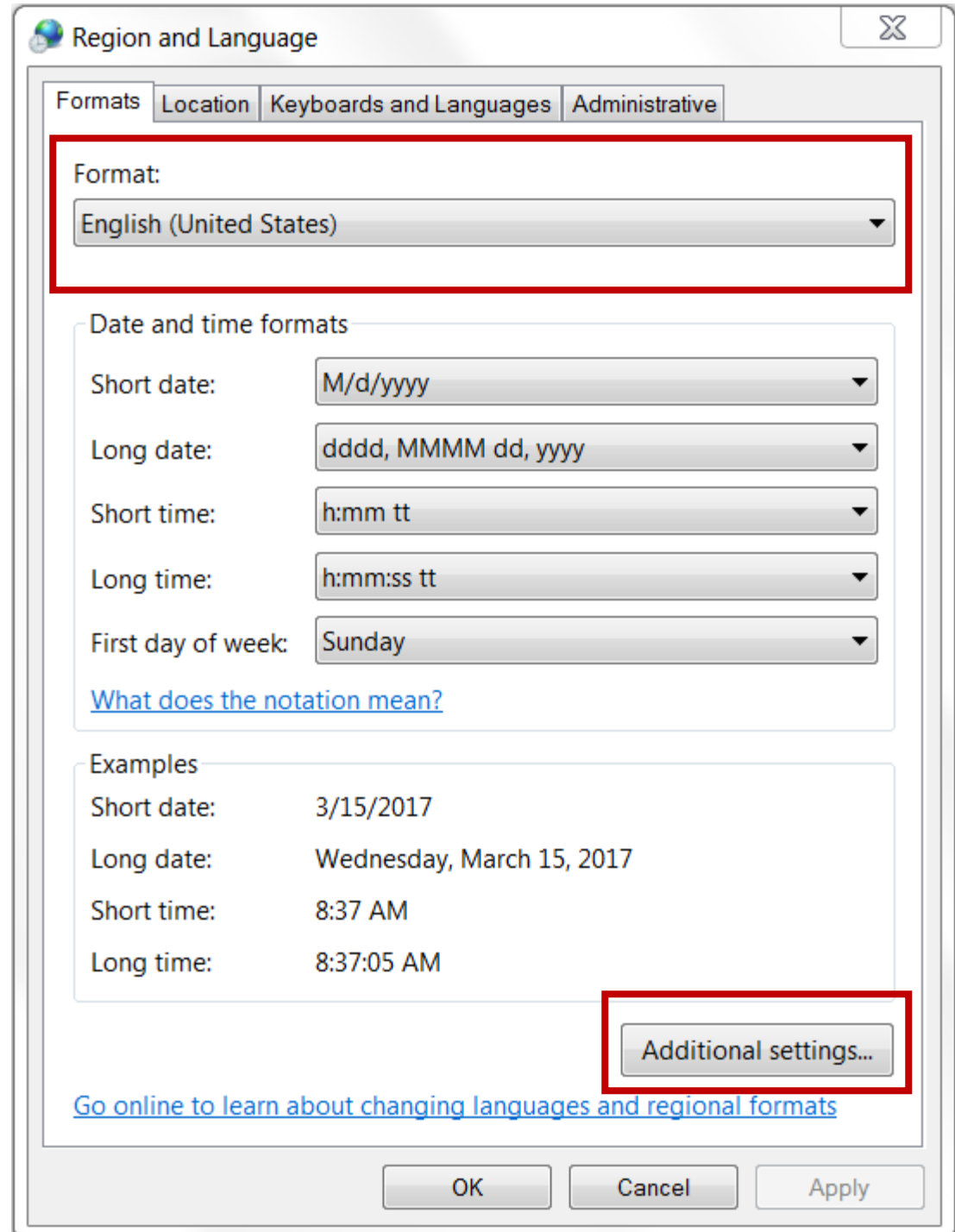
Standard digits: 0123456789

Use native digits: Never

Click Reset to restore the system default settings for numbers, currency, time, and date. Reset

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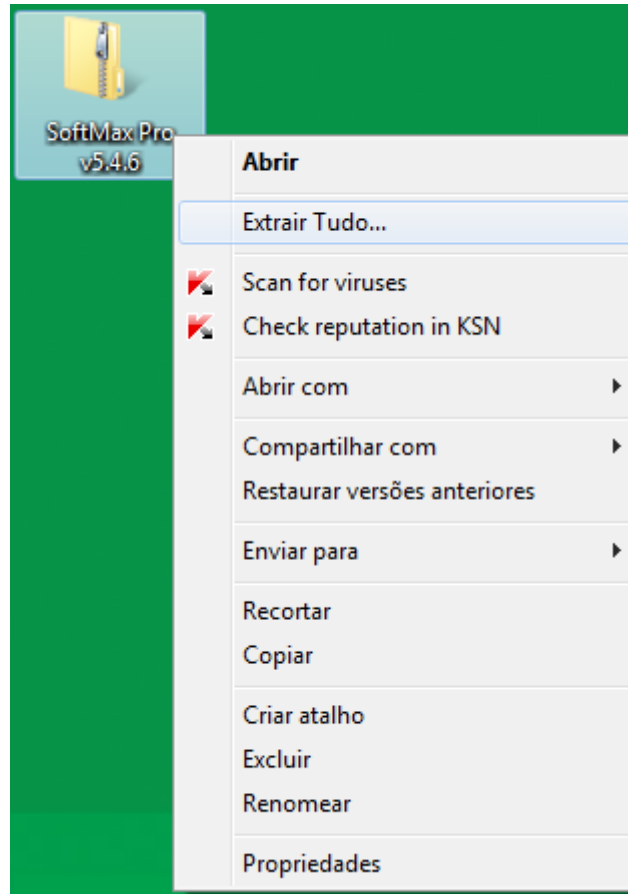
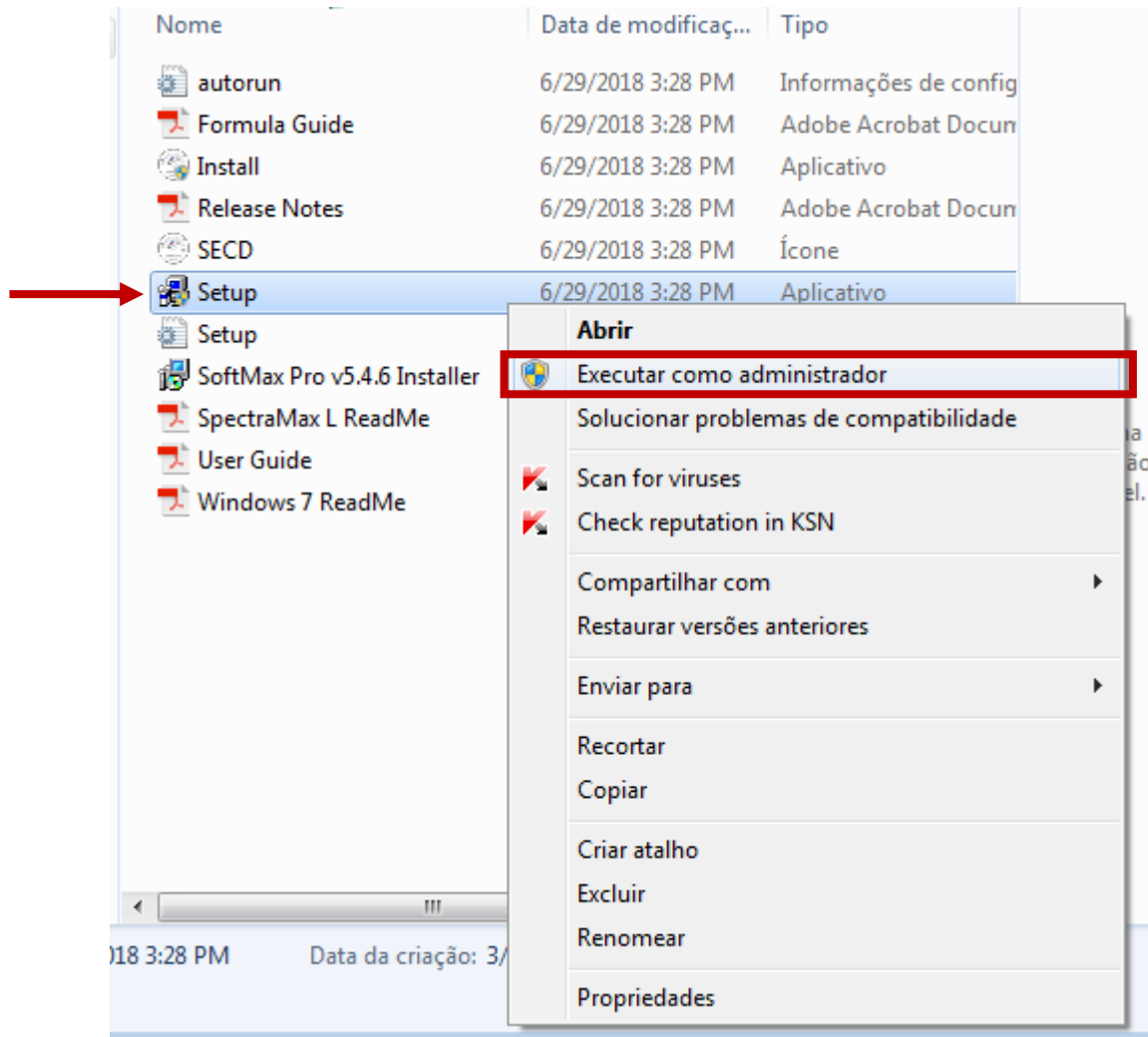
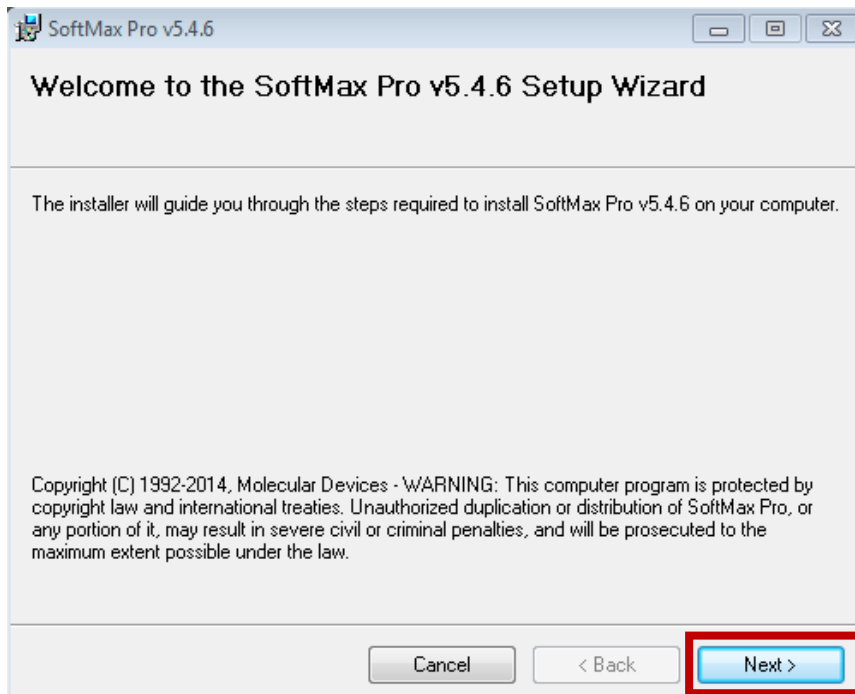
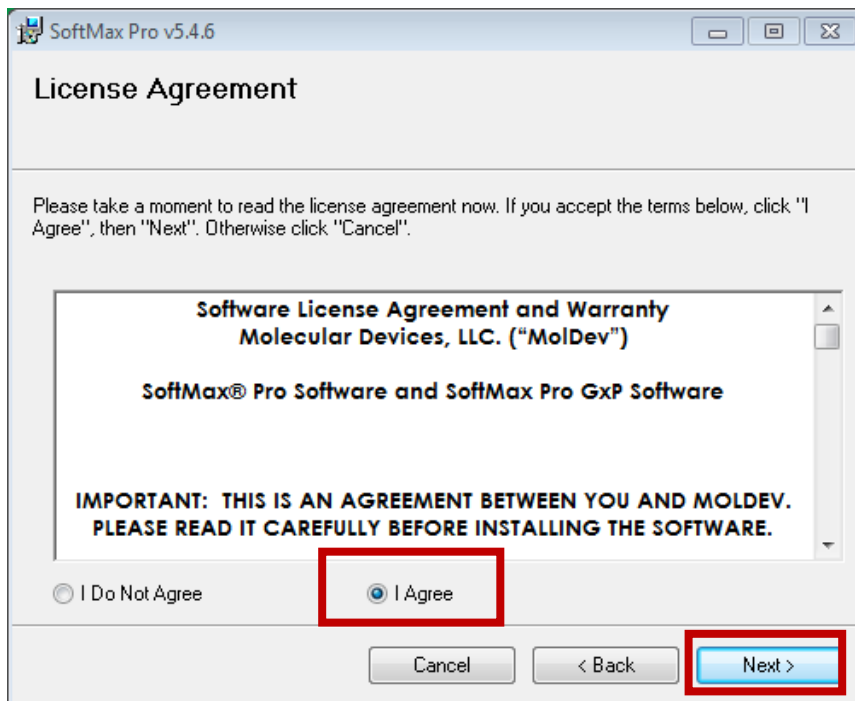
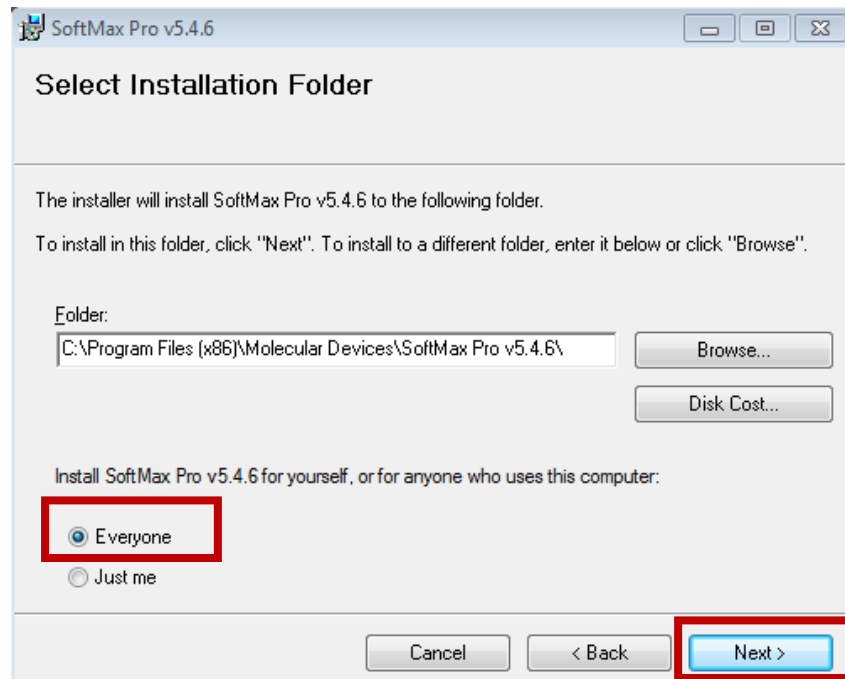
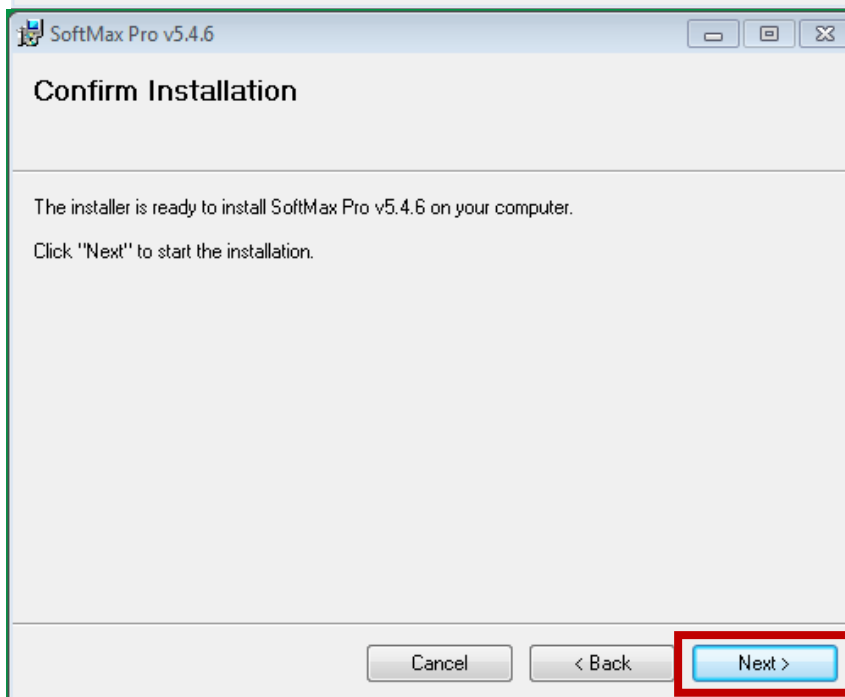
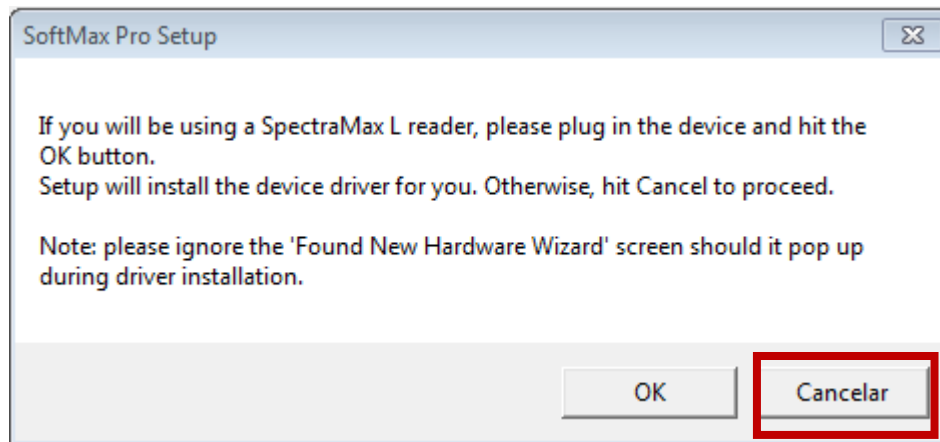
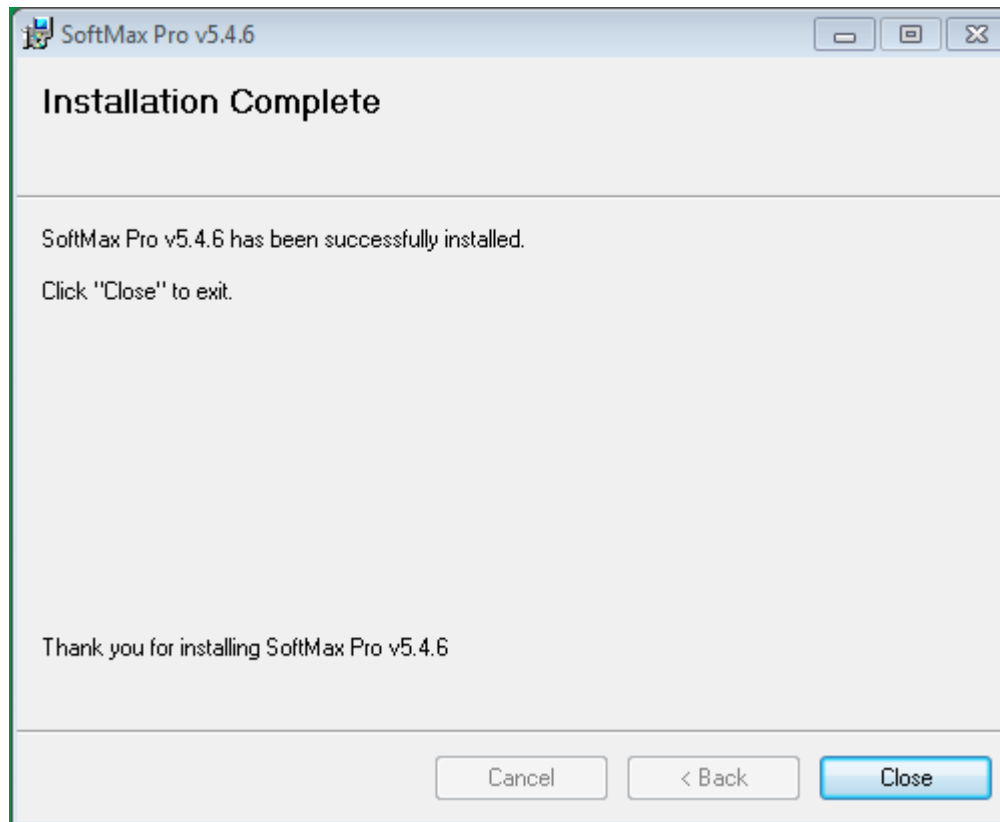
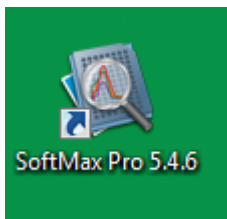
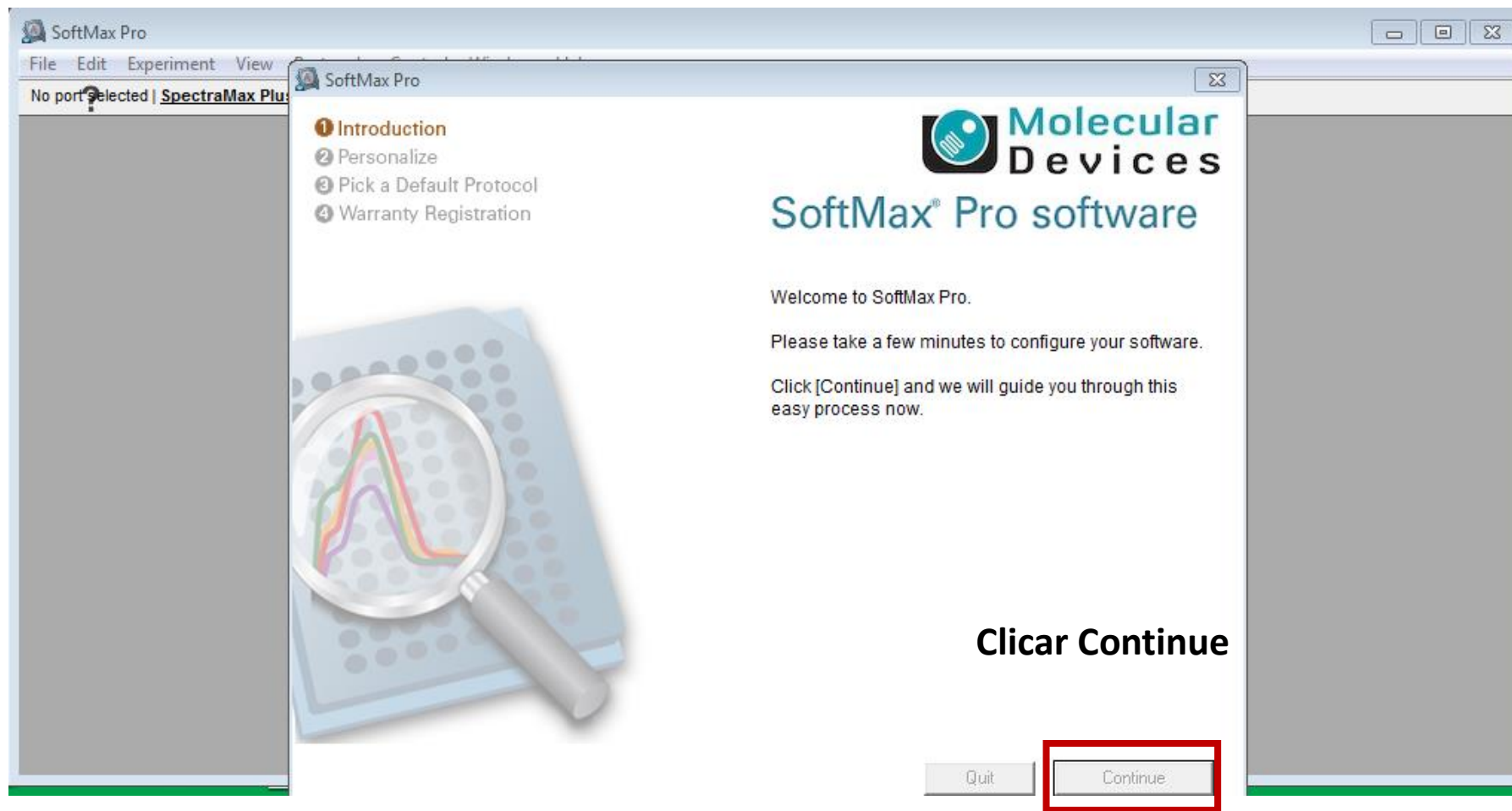
SoftMax[®] Pro software

You are required to personalize your copy of SoftMax Pro.

This required information will be used to identify your copy of the software.

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SoftMax[®] Pro

Molecular Devices

Before using SoftMax Pro software, please enter the following required information:

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The screenshot displays the SoftMax Pro software interface. The top menu bar includes File, Edit, Experiment, View, Protocols, Control, Window, and Help. Below the menu bar, a status bar shows 'No port selected' (highlighted with a red box), 'SpectraMax Plus', and 'StakMax' buttons. The main window is titled 'Untitled' and contains a tree view with 'Experiment#1', 'Notes#1', and 'Plate#1'. The 'Plate#1' section shows a 12x8 grid (columns 1-12, rows A-H) and a settings panel on the right. The settings panel includes 'Endpoint' (Lm1 450), 'Automix: Off', 'Calibrate: On', and 'Plate Last Read: No Data'. Below the grid, the text 'Wavelength Combination: !Lm1' and 'Data Type: Absorbance' is visible. The bottom of the window shows a 'CuvetteSet#1' section with 'Settings', 'Template', 'Reduction', and 'Display' buttons.

Plate#1												
	1	2	3	4	5	6	7	8	9	10	11	12
A												
B												
C												
D												
E												
F												
G												
H												

Endpoint
Lm1 450
Automix: Off
Calibrate: On
Plate Last Read:
No Data

Wavelength Combination: !Lm1
Data Type: Absorbance

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